




Stability –
giving you a reliable research tool
you can count on.

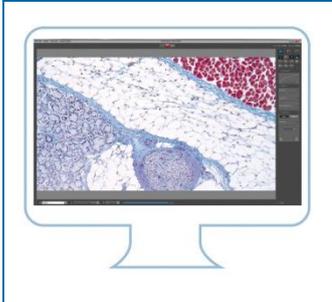

Transfer your software settings

In case you will transfer the settings of PROGRES GRYPHAX® software to another computer workstation you can copy the settings file from old computer to the new one.


Furthermore, in case of trouble shooting with the PROGRES GRYPHAX® software, you can rename the file and restart the software with default settings again.

To do so you need to copy or rename the settings file: "lastsession.ini" under:

Windows OS	"C:\ProgramDataJenoptik\PROGRES GRYPHAX\profiles"
Mac OS	"MacHD\Library\Application SupportJenoptik\PROGRES GRYPHAX\..."
Linux OS	"/var/lib/Jenoptik/PROGRES GRYPHAX/profiles/..."


For WINDOWS operation systems please follow the steps below to copy or rename PROGRES GRYPHAX® software settings file:

(1) Go to the control panel and activate: "show hidden folders" on your computer

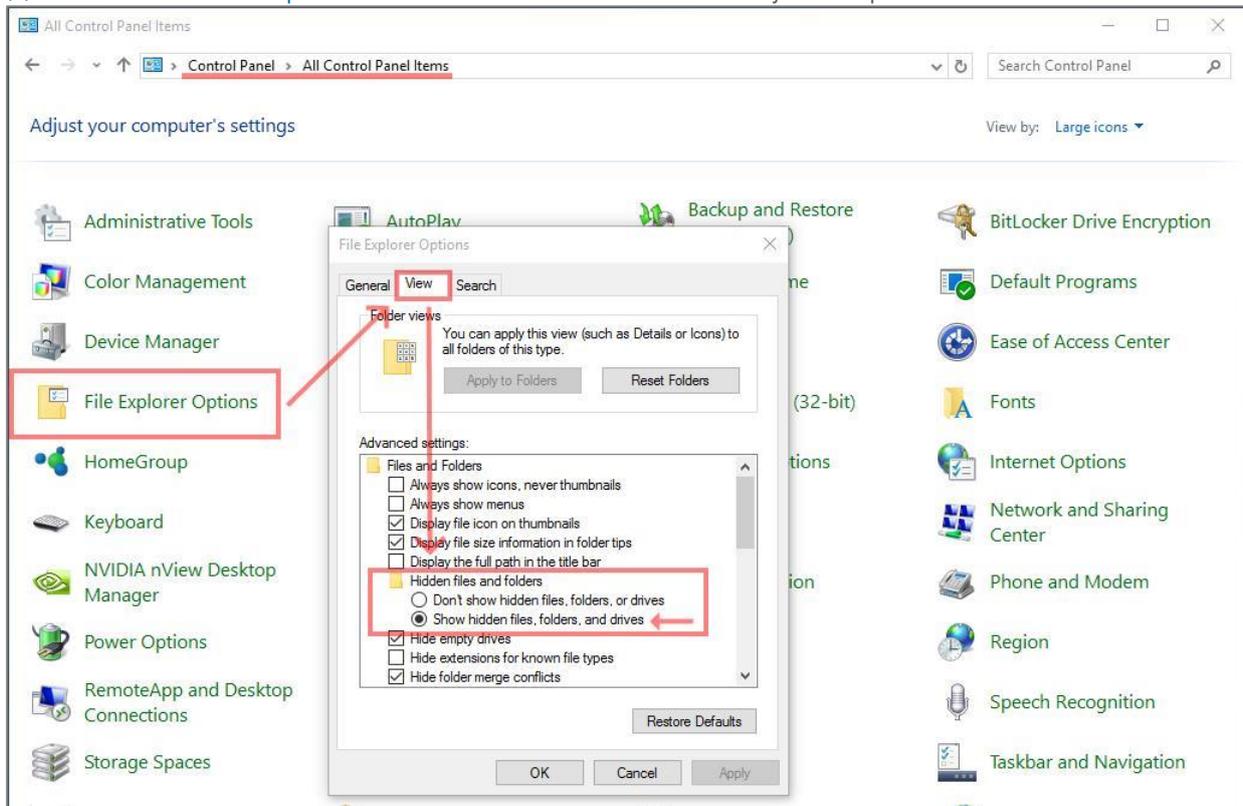




(2) Navigate to drive: "C:" and search for hidden folder: "ProgramData"

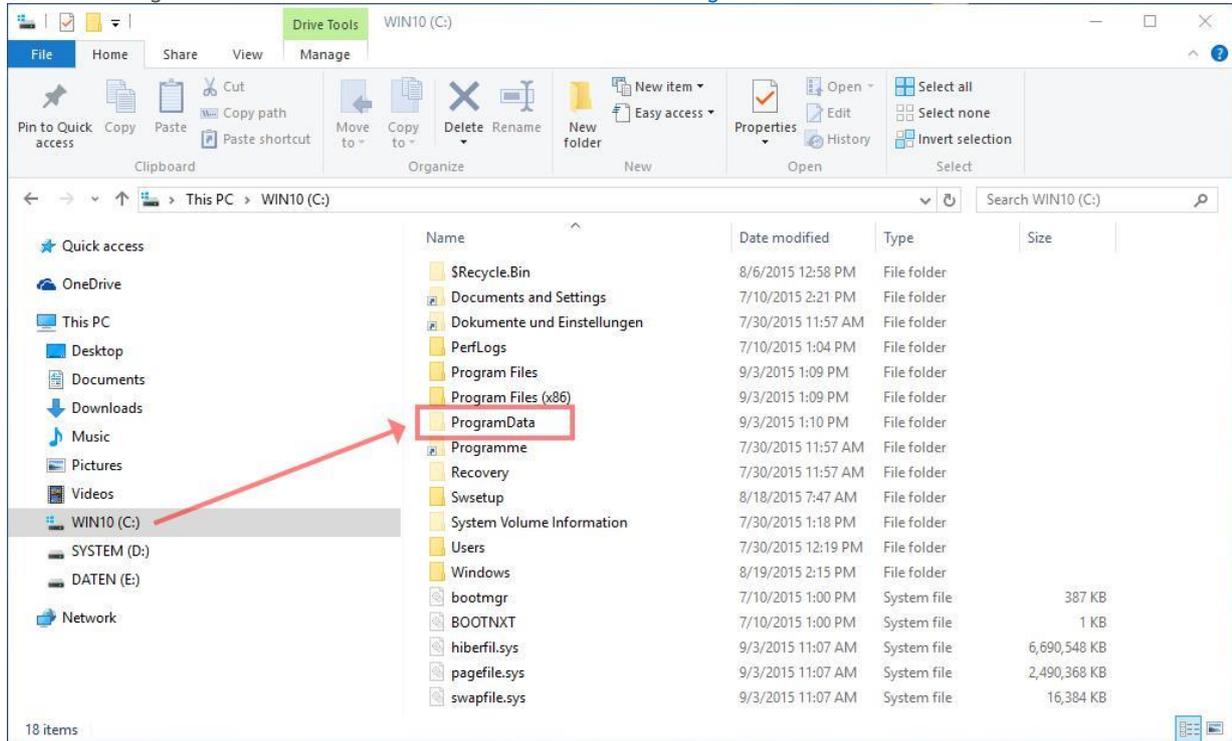

(3) Rename or copy settings file: "lastsession.ini" under: "C:\ProgramData\Venoptik\PROGRES GRYPHAX\profiles"

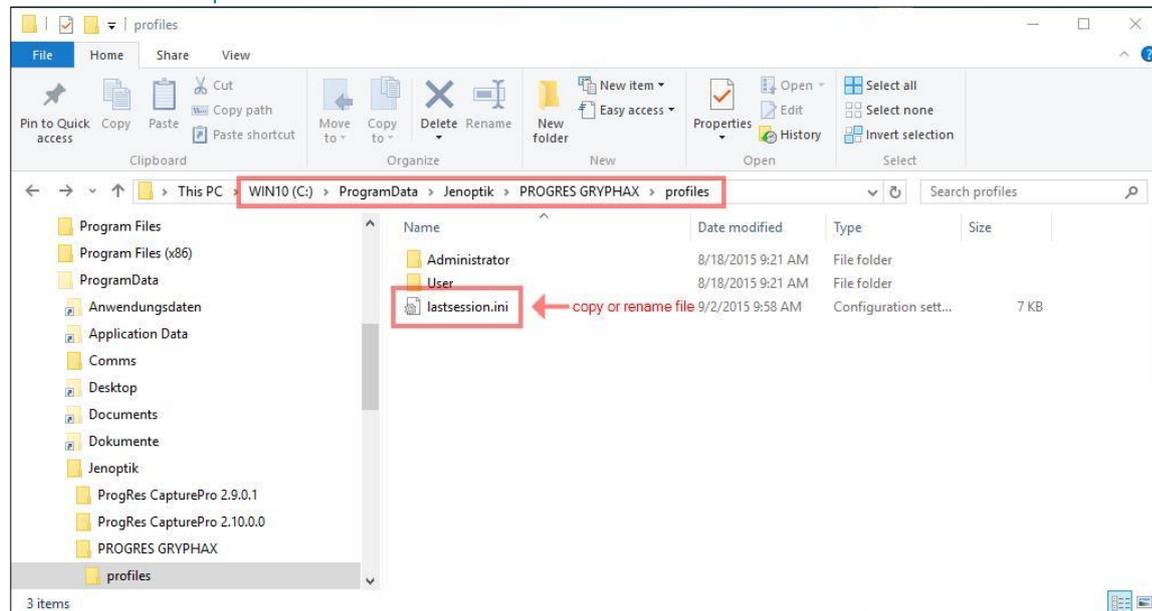

Please note again: Delete or rename of settings file "lastsession.ini" will **lose all** previous software settings and default settings will be used instead!

Your PROGRES GRYPHAX® Team.



JENOPTIK | Healthcare & Industry
JENOPTIK Optical Systems GmbH
Göschwitzer Straße 25