




### 33. Keyboard shortcuts

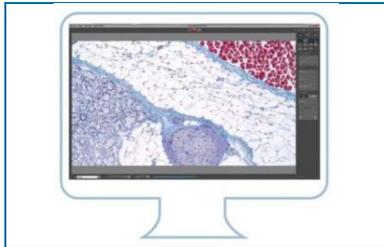

## Keyboard Shortcuts of JENOPTIK GRYPHAX<sup>®</sup> software

JENOPTIK GRYPHAX software is **workflow-optimized** and very easy to use. Therefore, software includes many **time-saving** keyboard shortcuts.

Keyboard button:	Shortcuts for:
F1	Open <b>help</b>
F2   ENTER	Start <b>capturing</b> in "Capture Mode" (single shot, time-lapse, video)
F3	Return to " <b>Live</b> " mode
F4	Open <b>Individual Save</b> dialog
F5	Activates   deactivates live image <b>freeze</b>
F6*	Switch between <b>Device configurations</b> at Status bar
F7*	Switch between <b>Magnifications / Lens</b> calibrations at Status bar
F11   F	Switch between " <b>Presentation view</b> " and "Standard view"
DEL	<b>Delete</b> selected images from "Gallery" or marked drawings (measurement   annotation)
Z	Switch <b>zoom</b> between "1:1" mode and "Fit to screen" mode
CTRL** + P	<b>Print</b> selected image(s) from "Gallery"
CTRL** + A	<b>Select all</b> (all thumbnails from "Gallery")
CTRL** + G	Open   close " <b>Gallery</b> "
CTRL** + O	Open " <b>Storage Options</b> " to change destination folders for gallery (images, videos)
CTRL** + T	Open   close " <b>Toolbar</b> "
CTRL** + D	Open   close " <b>Treeview</b> " to change and manage media destination folder fast & easy
G   C*	Activates   deactivates " <b>Grid</b> " or " <b>Crosshair</b> " view
V*	Activates   deactivates Image comparison " <b>Split Image view</b> "
ESC	<b>Leave</b> "Presentation view"   <b>remove</b> drawings during creation (measurement   annotation)
Arrow up   down	Gallery <b>thumbnail</b> up   down
Page up   down	Gallery <b>page</b> up   down
Home	<b>Jump</b> to "Gallery" <b>first</b> page
End	<b>Jump</b> to "Gallery" <b>last</b> page
Insert   Ins*	to insert user defined layer image while <b>Z-Stack   EDF</b> image record is running
SPACE BAR	Activates " <b>Magnifier</b> " or " <b>Hand tool</b> "
+   -	to <b>zoom IN</b> or <b>OUT</b> while "Magnifier" is active

\*valid as of version 2.3.0 | \*\*for Mac users: CMD instead

**Hint:** We all want to boost our productivity, so we recommend printing out or saving this list of keyboard shortcuts to enhance user's daily work.